

Dilucup Elegance BPW (Buffered Peptone Water)

Use and description: Ready-to-use plastic cups, each contains 9 mL medium to be used for diluting microbiological samples together with a special shaker, Dilushaker or Diligent Shaker. By adding a sample of 1 mL to the first cup a dilution of 1:10 will be achieved. 1 mL from this cup is then transferred to the next to get a dilution of 1:100. The procedure is continued until the required dilution is reached.

Composition per litre:

Peptones	10,00 g
NaCl	5,00 g
Buffers	3,56 g

pH: 7,0 +/-0,2 at 25°C

Volume per cup: 9 mL +/- 2 %

Sterilization: E-Beam

Material of the cups: PET

Expiry date: When unopened 30 months after production.

Storage: Unopened: +2 to +30°C. +2 to +8°C when the box is open.


Packaging: The cups are delivered in blisters of 7 pre-cut rows of 3 cups (40DEBPW3) or 6 cups (40DEBPW6). Each box contains 672 cups distributed in 2 bags of 336 cups. Weight and volume: 8,1 kg/box - 0,032 m³/box

DILUCUP ELEGANCE BPW 2 X 16 X 21 CUPS

ARTICLE NUMBER: 40DEBPW3

DILUCUP ELEGANCE BPW 2 X 8 X 42 CUPS

ARTICLE NUMBER: 40DEBPW6

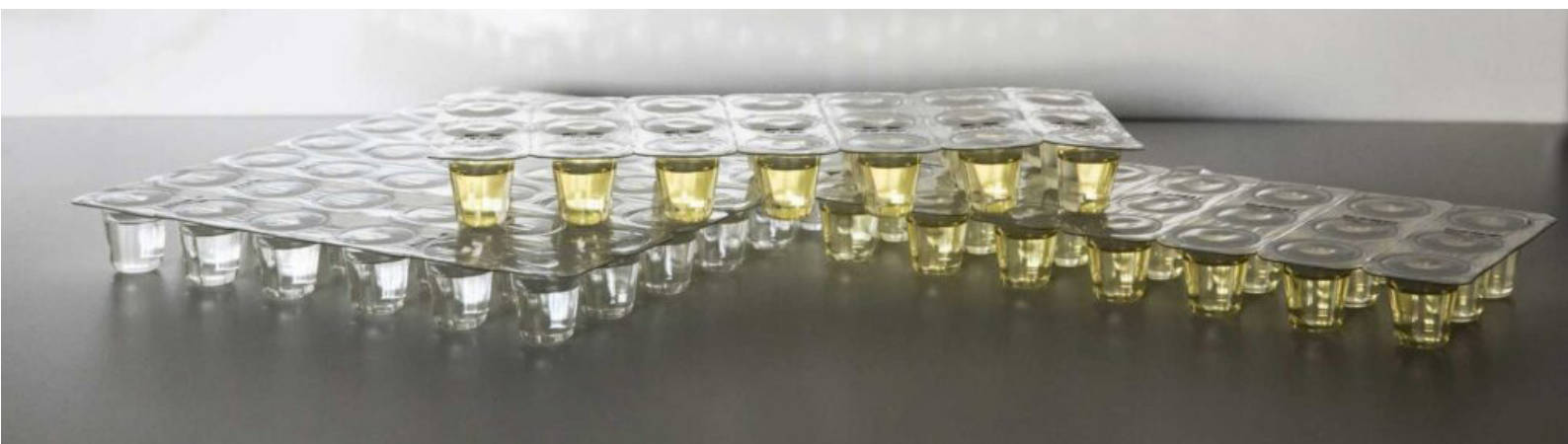

Dilucup Elegance BPW (Eau Peptonée Tamponnée)

Description & utilisation: Cupules de polyéthylène prêt-à-l'emploi contenant 9 mL de diluant. A utiliser sur un agitateur dédié (Dilushaker ou Diligent Shaker) pour la réalisation de dilutions décimales en microbiologie alimentaire. L'ajout d'1 mL d'échantillon dans une première cupule permet la réalisation d'une dilution 1:10. Le transfert d'1 mL de cette première cupule dans la cupule suivante permet la réalisation d'une dilution 1:100. Cette procédure doit être répétée jusqu'à l'obtention du niveau final de dilution souhaité.

Composition (par litre):

Peptones	10,00 g
NaCl	5,00 g
Tampons	3,56 g

pH: 7,0 +/-0,2 à 25 °C

Volume par cupule: 9 mL +/- 2 %

Stérilisation: Par E-Beam (faisceau d'électrons)

Matériau: PET

Date d'expiration: 30 mois à partir de la date de production.

Conservation: Entre +2 et +30°C avant ouverture. Entre +2 et +8°C ensuite.

Conditionnement: Les cupules sont présentées en blisters de 7 rangées pré-découpées de 3 cupules (40DEBPW3) ou 6 cupules (40DEBPW6). Chaque boîte contient 672 cupules réparties en 2 sachets de 336 cupules.

Poids et volume: 8,1 kg/boîte - 0,032 m³/boîte

DILUCUP ELEGANCE BPW 2 X 16 X 21 CUPULES

CODE ARTICLE: 40DEBPW3

DILUCUP ELEGANCE BPW 2 X 8 X 42 CUPULES

CODE ARTICLE: 40DEBPW6

