

interscience



Anabac[®]

Autoclave deodorant

Give some air to your lab!

Laboratories are very concerned about hygiene cleanliness and comfort of work. Try **Anabac®**, autoclave deodorant, for a perfumed touch in your lab. **Anabac®** autoclave deodorant bring the discrete final touch, the refinement of cleanliness.



Classic

Well perfumed
All applications

- *mint and eucalyptus fragrance*



Poma

Sweet perfume
Food industry

- *apple fragrance*



Citrus

Sweet perfume
Food industry

- *lemon fragrance*



Floral

Well perfumed
All applications

- *flower fragrance*




Peach

Well perfumed
All applications

- *peach fragrance*




1.

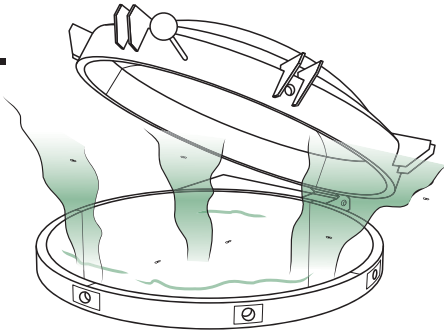


During sample sterilization, bad smell can emerge from the autoclave.

2.

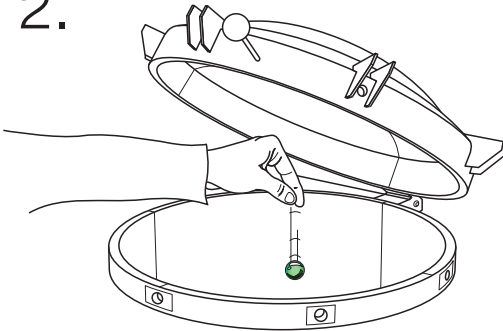

Place one capsule (or 2 if the odor is strong) in the water at the bottom of the autoclave.

3.

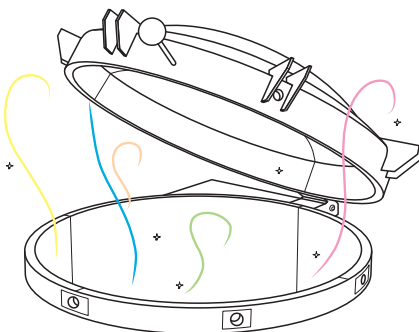

After a sterilization cycle, the capsule melts, allowing the perfume to be released.

Advantages

- *Work comfort*
- *Compatible with any autoclave*
- *No effect on the autoclaving*
- *Pleasant perfume*
- *100% biodegradable*

Applications

- For any kind of autoclave (max: 200 L), recommended for food industry.

Specifications

- Pot of 100 capsules
- Pot weight: 160 g
- 5 fragrances: - Classic (ref: 320 100)
- Poma (ref: 320 200)
- Citrus (ref: 320 300)
- Floral (ref: 320 400)
- Peach (ref: 320 500)



Please send this request to a local distributor or to interscience.

Anabac® autoclave deodorant

- Quotation
- Free "5 fragrances" sample

I want to receive informations on:

- BagMixer® 400 ml blender
- JumboMix® 3500 ml blender
- BagFilter® filter-bags
- BagTools® accessories
- Gravimat® dilutor
- easySpiral® Spiral plater
- Scan® 500 & 1200 colony counters
- Scan® 100 manual colony counter

Name.....

Company.....

Address.....

Zipcode/city..... Country.....

Phone..... Fax.....

Email.....

Web.....

interscience

INTERNATIONAL

30, Ch. Bois des Arpents - 78860 St Nom - FRANCE
T: +33 (0)1 34 62 62 61 - F: +33 (0)1 34 62 43 03
info@interscience.fr - www.interscience.fr

USA & CANADA

199 Weymouth ST - ROCKLAND - MA 02370 - USA
T: +1 781 792 2133 - F: +1 781 792 2134
info@intersciencelab.com www.intersciencelab.com

In addition



Clinet® 101

Chlorhexidine digluconate disinfecting wipes

Ideal for laboratories, hospitals, samplers, home use.



Clinet® 202

Individually packed chlorhexidine digluconate disinfecting wipes

Ideal for laboratories, hospitals, samplers, transports, hand hygiene.



Clinet® 303

Bactericidal and fungicidal spray

Ideal for large surfaces, hard to reach places, angle of walls, rough edges.



Clinet® 404 SaniFood®

Food contact approved disinfecting wipes

For food professionals: disinfection without rinsing of surfaces and materials in contact with food.



Clinet® Anti-algae

Algicide for laboratories water baths.

1 liter bottle with incorporated doser, Clinet®Anti-algae is used diluted at 0.1% to inhibit the growth of algae in water baths.

Anabac_B_0314 - Pictures and information in this leaflet are not contractually binding. Clinet® 101, Clinet® 202, Clinet® 303, Clinet® 404 SaniFood, Clinet® Anti-algae are registered trademarks.

