

# Dilucup Elegance NACL

**Use and description:** Ready-to-use plastic cups, each contains 9 mL medium to be used for diluting microbiological samples together with a special shaker, Dilushaker or Diligent Shaker. By adding a sample of 1 mL to the first cup a dilution of 1:10 will be achieved. 1 mL from this cup is then transferred to the next to get a dilution of 1:100. The procedure is continued until the required dilution is reached

**Composition per litre:**

NaCl 8,5 g  
Buffers 1,4 g

**pH:** 7,0 +/-0,2 at 25°C

**Volume per cup:** 9 mL +/- 2 %

**Sterilization:** E-Beam

**Material of the cups:** PET

**Expiry date:** When unopened 30 months after production.

**Storage:** Unopened: +2 to +30°C. +2 to +8°C when the box is open.


**Packaging:** The cups are delivered in blisters of 7 pre-cut rows of 3 cups (40DENACL3) or 6 cups (40DENACL6). Each box contains 672 cups distributed in 2 bags of 336 cups. Weight and volume: 8,1 kg/box - 0,032 m<sup>3</sup>/box.

DILUCUP ELEGANCE NACL 2 X 16 X 21 CUPS

**ARTICLE NUMBER:** 40DENACL3

DILUCUP ELEGANCE NACL 2 X 8 X 42 CUPS

**ARTICLE NUMBER:** 40DENACL6

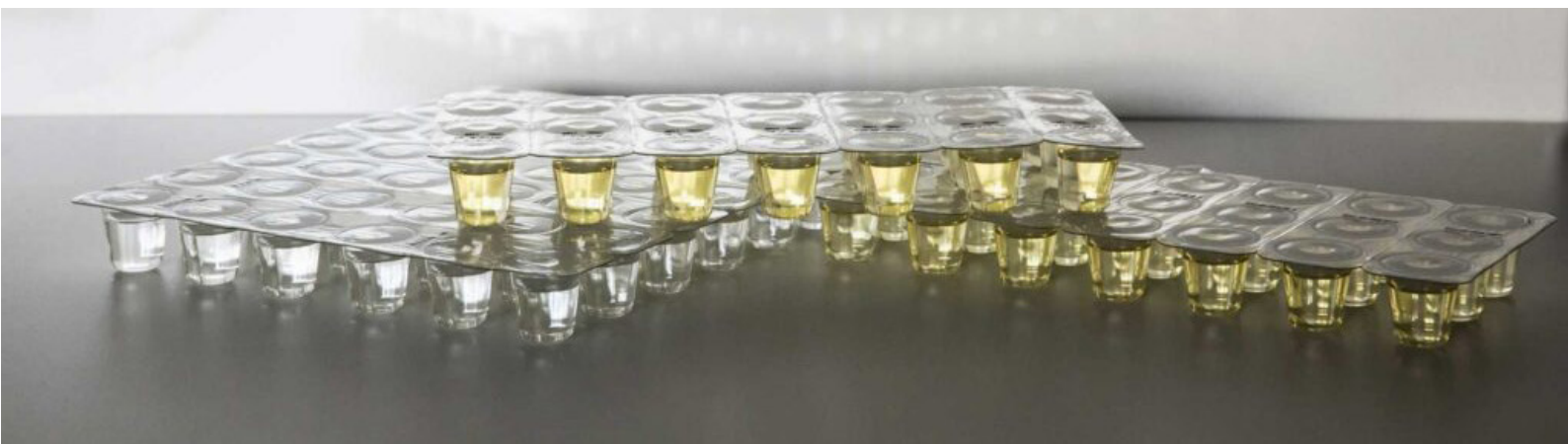

# Dilucup Elegance NACL

**Description & utilisation:** Cupules de polyéthylène prêt-à-l'emploi contenant 9 mL de diluant. A utiliser sur un agitateur dédié (Dilushaker ou Diligent Shaker) pour la réalisation de dilutions décimales en microbiologie alimentaire. L'ajout d'1 mL d'échantillon dans une première cupule permet la réalisation d'une dilution 1:10. Le transfert d'1 mL de cette première cupule dans la cupule suivante permet la réalisation d'une dilution 1:100. Cette procédure doit être répétée jusqu'à l'obtention du niveau final de dilution souhaité.

**Composition (par litre):**

NaCl 8,5 g  
Tampons 1,4 g

**pH:** 7,0 +/-0,2 à 25 °C

**Volume par cupule:** 9 mL +/- 2 %

**Stérilisation:** Par E-Beam (faisceau d'électrons)

**Matériau:** PET

**Date d'expiration:** 30 mois à partir de la date de production.

**Conservation:** Entre +2 et +30°C avant ouverture. Entre +2 et +8°C ensuite.

**Conditionnement:** Les cupules sont présentées en blisters de 7 rangées pré-découpées de 3 cupules (40DENACL3) ou 6 cupules (40DENACL6). Chaque boîte contient 672 cupules réparties en 2 sachets de 336 cupules

**Poids et volume:** 8,1 kg/boîte - 0,032 m<sup>3</sup>/boîte

DILUCUP ELEGANCE NACL 2 X 16 X 21 CUPULES

**CODE ARTICLE:** 40DENACL3

DILUCUP ELEGANCE NACL 2 X 8 X 42 CUPULES

**CODE ARTICLE:** 40DENACL6

