**U.S. English**

*Product Number:*  
8565

***Reveal<sup>®</sup> 3-D***  
*for Coconut*

# Reveal<sup>®</sup> 3-D for Coconut

Product Number: 8565

## Introduction

Reveal<sup>®</sup> 3-D for Coconut allergen is uniquely designed with three lines of detection and can be used virtually anywhere to screen environmental swabs, clean-in-place (CIP) rinses, food products and, ingredients for the presence of coconut. Its 3-D technology ensures greater reliability with screening than ever before.

## Intended Use

Reveal 3-D for Coconut allergen is intended for use only in an industrial food manufacturing/preparation context. Because of the problems of adequate sampling and extraction of coconut, it is not suitable for the testing of foods to be consumed in the home or in a restaurant by allergic individuals. The test detects coconut residue in CIP rinses, environmental swabs, and foods.

## Detection limit

The test uses highly specific antibodies to detect coconut. Coconut residues can be detected from various surfaces using the provided environmental swabs or from rinses and foods.

Utilizing the environmental swabs supplied, levels of 1.0 µg/100 cm<sup>2</sup> of coconut protein on surfaces can be detected.

When analyzing rinses or foods, coconut residues are detectable as low as 1 ppm coconut protein for rinses and 5 ppm coconut protein for foods. The presence of cleaners and sanitizers in rinses, as well the type of food matrix being tested, can affect limit of detection.

## Test Performance

Reveal 3-D for Coconut allergen has undergone rigorous validation to evaluate the specificity, sensitivity, robustness and intra- and inter-batch variability of the test method on CIP rinses, environmental swabs, and foods.

## Sample Compatibility

Reveal 3-D for Coconut allergen is designed to detect coconut on environmental surfaces and in CIP rinses, and foods. Although every effort has been made to validate as many variables as possible, there are some sample types that are not suitable for testing.

User should perform in-house matrix-specific spike and recovery validation work in conjunction with a validated laboratory assay (e.g. Veratox<sup>®</sup>) to help confirm Reveal 3-D for Coconut results. This process will highlight any problematic matrices encountered.

For certain complex food matrices, it is recommended that the use of a more broadly inclusive test method, such as Veratox<sup>®</sup> for Coconut allergen, be used. Existing commodity validations should be revalidated when suppliers or the manufacturing process has been changed. Please contact a NEOGEN representative or distributor for additional details.

## Materials Provided

Reveal 3-D for Coconut allergen (NEOGEN item 8565) includes:

1. Instructions for use
2. 1 foil pouch, containing 10 Reveal 3-D for Coconut allergen devices
3. 10 sachets containing 4 mL Type 8 Extraction Buffer
4. 10 sample tubes and caps
5. 10 individually packaged sterile swabs with break-off tips

## Materials Required for Food Extraction but Not Provided

1. 3-D Food Buffer (NEOGEN item 8504)
2. 50 mL graduated tubes for extraction (NEOGEN item 9381)

## Sampling Technique

1. CIP rinses and foods

As only a small amount of material is required for Reveal 3-D for Coconut allergen, it is important to test a representative portion of the sample.

2. Environmental swabbing

The swabs supplied are intended to be used for the collection of environmental samples from which the presence of coconut residues can be tested. This method can be used to validate the adequacy of cleaning and/or to identify problem areas (e.g., unwanted buildup of coconut residue in processing equipment).

## Precautions

1. For environmental testing in an industrial food manufacturing/preparation, or labeling enforcement context only.
2. Do not use any part of the test beyond the expiry date.
3. Do not open the foil pouch until just before use.
4. Ensure the foil pouch is sealed tightly after removal of device.
5. Always store the kit between 2–8°C (35–46°F). Avoid freezing.
6. Bring kit to room temperature 18–30°C (64–86°F) prior to use.

## Test Storage / Sample Stability

Store the Reveal 3-D for Coconut allergen kit between 2–8°C (35–46°F), and use within the expiry date stated on the outer label. Extracted samples should be tested within 3 hours of extraction.

## Limitations


A negative test from a surface or sample cannot exclude the possibility that the food contains coconut residue since it may be distributed unevenly on the surface or within the food and may be below the limit of detection of the test.

Reveal 3-D for Coconut allergen is qualitative and should only be used as a preliminary screen for the presence of coconut residue. The validity of results obtained should be viewed in conjunction with data from a quantitative assay, such as Veratox for Coconut allergen.


Hydrolyzed and fermented proteins may not be detectable with antibody-based test methods for allergen testing. Although the proteins may be undetectable in the assay because of their properties, there still could be allergenic protein residue present.

Validate detection by testing a positive control from the source of allergen or ingredient that is available in your manufacturing environment and that may pose an allergen contamination risk to ensure the test can detect the allergen of concern.

Contact NEOGEN Technical Services for additional information.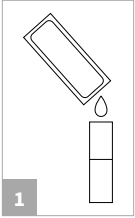

1

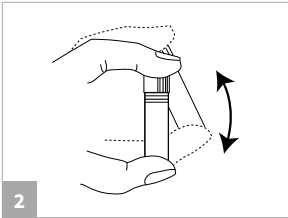
2


## Liquid Sampling

Remove the following and allow to equilibrate at room temperature before use (20–30 minutes out of refrigerator):

1. 1 sample tube and cap
  2. 1 Type 8 extraction buffer sachet
  3. 1 Reveal 3-D Coconut test device (in foil pouch)
1. Carefully tear/cut off the top of the Type 8 extraction buffer sachet and add the entire contents to the sample tube.
  2. Add 0.25 mL (250 ul) of sample to the sample tube.
  3. Secure the white cap and shake for 1 minute.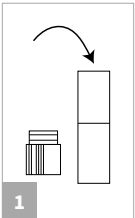

1

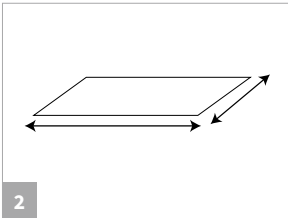
2

## Swab Sampling


For each sample to be tested, remove the following from refrigeration and allow to equilibrate at room temperature before use (20–30 minutes out of refrigerator):

- 1 sample tube
  - 1 Type 8 extraction buffer sachet
  - 1 Reveal 3-D for Coconut allergen device (in foil pouch)
  - 1 sterile swab
1. Carefully tear/cut off the top of the Type 8 Extraction Buffer sachet and add the entire contents to the sample tube.
  2. Estimate a swabbing area of approximately 10 x 10 cm. Alternatively, use the swab to collect samples of contamination from problem areas (e.g., processing equipment, filler heads).
  3. Gather the sample with the swab, using one of the following methods: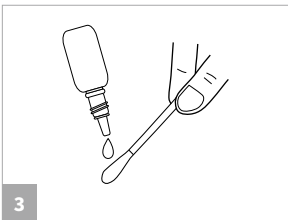

3


For dry surfaces: Remove a sterile swab from the packaging and wet with extraction buffer in the sample tube. Swab a 10 x 10 cm area by using a crosshatch technique revolving the swab on the surface. Repeat this swabbing procedure using movements at right angles to those used in the first swabbing.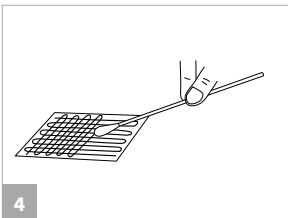


4

For wet surfaces: Remove a sterile swab from the packaging and swab a 10 x 10 cm area by using a crosshatch technique revolving the swab on the surface. Repeat this swabbing procedure using movements at right angles to those used in the first swabbing. Do not moisten swab prior to use.


4. Return the swab to the extraction buffer in the sample tube and carefully break off the moistened end at the pre-scored mark so that it remains in the tube.
5. Secure the cap of the sample tube, taking care to ensure the stem does not prevent the tube from being properly sealed. Shake for 1 minute.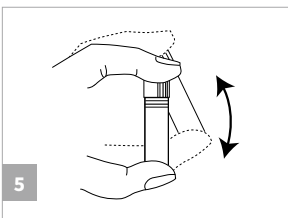

5

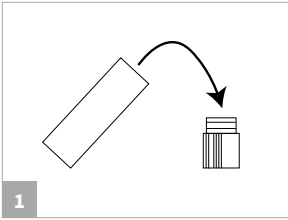1

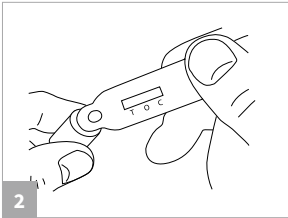
2


## Swab and Rinse Testing

1. Remove the lid and fill it with the liquid from the tube. Any froth should remain in the tube.
2. Dip the head of the Reveal 3-D device into the liquid in the lid. Ensure the cavity is saturated with the liquid.
3. Leave the cavity saturated until liquid is seen running in the test window.
4. Place device on a flat surface and allow test to develop for 5 minutes.


**Note:** It is essential to place the device flat on a level surface as soon as the liquid has entered the test window to stimulate flow through the device. Additionally, the devices are pre-stripped with pale green loading dye in positions T (test), O (overload), and C (control). The loading dye is removed from the test window as the sample flows through the device.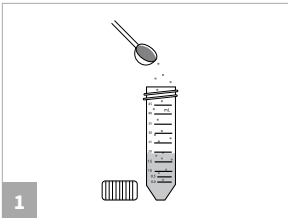

1

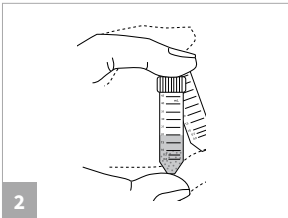2

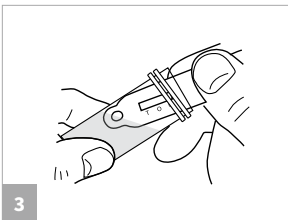3

## Food Extraction and Testing

The extraction of food samples requires the use of NEOGEN's 3-D Food Buffer (NEOGEN item 8504).





From the test kit, remove the appropriate number of the following and allow to equilibrate at room temperature before use (20–30 minutes out of refrigerator):

3-D Food Buffer

Reveal 3-D for Coconut device (in foil pouch)

1. Add 20 mL of 3-D Food Buffer into a new 50 mL tube (18 mL for liquid samples) or a similar container.
2. Weigh out 2 g (2 mL for liquid products) of sample and add to tube with buffer.
3. Secure the caps and invert to suspend sample. Shake vigorously or vortex for 1 minute.
4. Tilt the tube at approximately a 45 degree angle. Dip the head of the Reveal 3-D device into the liquid. Ensure the cavity is saturated with the liquid, and that no liquid gets on the test window.
5. Leave the cavity saturated until liquid is seen running in the test window
6. Place device on a flat surface and allow test to develop for 5 minutes.

**Note:** It is essential to place the device flat on a level surface as soon as the liquid has entered the test window to stimulate flow through the device. Additionally, the devices are pre-stripped with pale green loading dye in positions T (test), O (overload), and C (control). The loading dye is removed from the test window as the sample flows through the device.

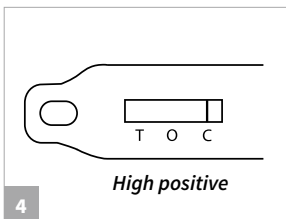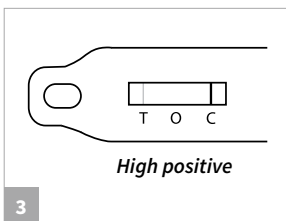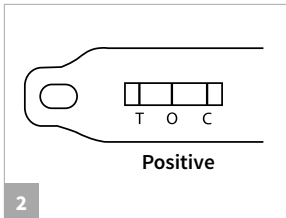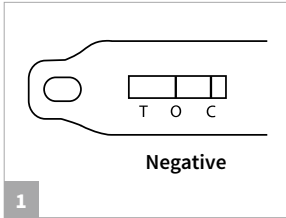## Reading Results

Read the result 5 minutes after dipping. Observations after 6 minutes may be inaccurate due to overdevelopment of the device.

1. Negative result  
No line at position T (test): Level of coconut below detection limit. (See limitations section).
2. Positive result  
Any intensity of line at position T (test): Level of coconut above detection limit.
3. High results  
No line is visible at position O (overload) and a line is faintly visible or absent at position T: Sample is overloaded with coconut.
4. Invalid results  
If no line appears at position C (control), then the test is invalid.

## Customer Service

NEOGEN Customer and Technical Services can be contacted through [NEOGEN.com](https://www.NEOGEN.com) and product training is available by request.

## SDS Information Available

Safety data sheets are available for all test kits at [NEOGEN.com](https://www.NEOGEN.com) or by calling 800.234.5333 or 517.372.9200.

## Terms and Conditions

NEOGEN's full terms and conditions are available [online](#).

## Warranty

NEOGEN makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, NEOGEN will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. NEOGEN shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

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